

Hans P. Harmon¹, Eric L. Akers¹, David A. Johnson¹, Erik Naylor¹, Jerome Clasadonte² & Philip G. Haydon² 1 - Pinnacle Technology Inc., Lawrence, KS 2 - Tufts University, Boston, MA

Introduction

- Because large variations in EEG waveforms are found between animals due to such parameters as screw placement, brain morphology and animal size a majority of researchers still rely on human visual analysis of EEG data for accurate interpretation of seizure events.
- Analysis of long EEG datasets is very resource intensive, creating a hindrance to large-scale projects.
- To help streamline data analysis, we have developed a software toolset for rapid analysis of EEG-based data.
- The software includes detailed power spectrum analysis with adjustable Hanning and Hamming windows along with customizable power band selection for all channels. Once power spectra have been calculated for a specific dataset, colorized heat maps and power frequency plots are available for visual analysis.
- Rapid identification of seizures within large datasets is accomplished based on parameters such as seizure length, specific frequency, root mean squared (RMS) power and line length.
- Identified seizure events can further be classified using embedded video and rated using standard Racine's Scale parameters. After identification has been completed, aggregated seizure events are analyzed according to number, severity, frequency, and circadian distribution over user-defined periods and adjustable for light/dark cycles.

Video Recording

- Video can be recorded at varying resolutions (320 x 240, 640 x 480), adjustable frame rates and from different angles to meet the needs of any experiment including chronic experiments lasting weeks (Figure 1).
- Video frame to data point accuracy is less than 100 ms.
- Individual seizure events can be isolated and exported with video recording as .avi video clips for use in presentations and papers.

Custom Plot Generation

• The seizure analysis tools provide a method of viewing and comparing various frequency band and line length data as well as the raw EEG signal on a single time plot. Various filters and scaling can be applied to the waveforms to help visualize the data (Figure 2).

Posit Guide	tion(y,t) e Locatio	03/30/2011 02:27:20 n 11/26/2012 05:04:11	5. 143 AM 🐳 🖬	20217			
F EMG	Full		Add a	20 <u>0</u>	the leave a		
Dat	ta Chann	els	f		/M NIM		
		EMG_Full	A 194		. I'L A DW MAA	AA A AMANA AMA	M. H. M
xis		🔽 Median Filter	11 pt	150	And man harmal "	MANNAN ULAN III	MANN W
6		💟 Mean Filter	7pt 🗧	- W	where a mode shift with	a data dha Andah Al	Contraction of the second second
xis		🔣 Bin	10 pt 6				
		Show Zero Line					
		Derivative	101	10 <u>0</u>	N 1991	۱ L	
		None					A. h
		1st Order	E		M man in	Warda (Vm)	my al
		2nd Order		50 V		Why I	Nov
		StdDev Bounds 2.0	00 🗧 🕻 🛯				
		🖻 EEG1_Theta				$\mathbf{\omega}$	U
		Median Filter	10 pt	۰			
		Mean Filter	6 pt				
		🔣 Bin	10 pt				
		Show Zero Line		-50	MARKER R. J. MARKER	A MALL AND A MALLAN	Minda Mari Jan L M
		Derivative			A MAL AM MAL AND A	an why have a like of	
		None					
		🔘 1st Order		-100		Y.	
		2nd Order	int				
		StdDev Bounds 2.0	00 🚔 [Ci.				
•			•	-152.4 03/2	29/2011 11:30:18.024 PM	103/30/2011 01:39:55 AM	103/30
		12: 		and the second se		Depo	1

A Software Toolset for Rapid Analysis of EEG Seizure and Video Data

▶ 🖩 🔀 🖓 🕫 🖉 👰 🙊 💯 🗘 📮 🛊 📮 🖉 🖉 👰 🙊 🖉 🖉 🖉 06/26/2011 01:44:21 PM

Figure 3: Seizure recording with integrated, synchronized video playback

- The seizure analysis software displays the data, annotations, and a synchronization line for video (Figure 4).
- Frequency data can also be displayed and is updated in real time based on the data signal
- A heat map of the data spectrum plot displays time seizure frequency Of for rapid red activity identification of seizure а event.



Seizure Analysis







Seizure and Video Synchronization



Figure 4: Raw EEG Data (top), Spectral Plot Data (bottom left), and Spectrum Heat Map Plot (bottom right).

- The seizure selection tool extracts database.
- The extracted features are then used to match patterns and identify similar seizures in other data sets (Figure 6).
- Features extracted include RMS power of a specific frequency band or line length.
- The features and the seizure waveform itself are stored in a database for future use in new analysis.

	e Search Setting	s Method	Method			
Siezure Type II	 Window Size 3.00 s 	Line Length	Line Length 👻			
View	Step Size 1.00 s	Eine Lei	Line Length Settings			
	Match 90.0 %	Threshold	Threshold 35605.70 len/sec Current Line Length		ec 主	
		Current Line				
1.7				-		
Search F	or Seizures			III Pa	use Search	
Current Locat	ion: 03/30/2011 08:59:5	57.000 AM				
Matched Sei	zures				🕕 Go To	
ID	Start Time	End Time	Value	Тур	Accept	
1 0	03/29/11 11:11:08	03/29/11 11:11:33	36589.7		Accept All	
2 1	03/30/11 02:21:31	03/30/11 02:21:51	38418.2		Remove	
3 2	03/30/11 03:05:52	03/30/11 03:06:16	41285.2		Merge Up	
		e:			' Merge Dowr	
					Next	
					Previous	
				5	Save TSV	
				01 1422		

a matched pattern in the EEG waveform (right).

- Seizure detection was performed using files. All files were compared to an expert's analysis (Figure 8).
- EEG and EMG seizure events.
- events within large-scale datasets.



Seizure Selection

features from an existing seizure and places the event and feature set into a



Figure 6: Seizure selected (in blue) and added to a pattern matching database. Frequency selection is shown in red.

Seizure Matching

- The seizure matching tool uses the seizure database to find similar seizure events and list them for human review (Figure 7).
- Pattern matching includes, but is not limited to, RMS power and line length.
- All the events can be accepted, rejected, or merged with nearby events to better classify seizures.



Figure 7: Pattern matching a seizure event. A list of matched events (left). A stored seizure event (middle). And

Testing

RMS power and line length analysis separately on five individual mouse data

AUGUKAUY UF SEIZUKE DETECTIU		ACCU	IRACY	OF	SEIZURE	DETECTIO	
------------------------------	--	------	-------	----	---------	----------	--

	# of Seizure Events Marked	Accuracy vs Expert Scorer	# of False Positives
Expert Scorer	21	N/A	N/A
RMS Power	23	100%	2
Line Length	21	100%	0

Figure 8: Seizure detection compared to manual scoring

Conclusions

• This software package provides comprehensive toolset for analysis and evaluation of

• Tools for automated seizure detection can be used for rapid identification of seizure

• Automated summary analysis tools provide researchers with rapid assessments of their data in both graphical and ASCII formatted outputs.